

TWO STEPS FOR FAST, ACCURATE PLATE READING



CLASSIFICATION ProtoCOL 2 Batch creation

- Differentiate between colour, size and shape
- Upper and lower count limits can be set i.e. <10cfu or >300cfu
- Selection of a counting area i.e. whole filter or half filter
- Separation of touching colonies
- Exclusion of unwanted items such as product
- Grid classification

MEASURE Count using ProtoCOL 2

- Automated count in seconds
- Detection of organisms as small as 43µm
- Average multiple plate counts
- Counts and images stored automatically
- Manually, add or delete colonies with an audit trail to comply with GMP/GLP
- Results can be directly transferred to a LIMS system, Excel or entered into one of ProtoCOL 2 customisable reports



A DIVISION OF THE SYNOPTICS GROUP


Synbiosis USA Headquarters:
5108 Pegasus Court Suite M
Frederick MD 21704 USA
Tel: 800-686-4451/301-662-2863
Fax: 301-631-3977
email: ussales@synbiosis.com

Synbiosis Europe and
International Headquarters:
Beacon House Nuffield Road
Cambridge CB4 1TF UK
Tel: +44 (0)1223 727125
Fax: +44 (0)1223 727101
email: sales@synbiosis.com

Website: www.synbiosis.com


A.022.11.10 All trademarks acknowledged

MICROBIAL LIMITS TEST - MEMBRANE FILTRATION

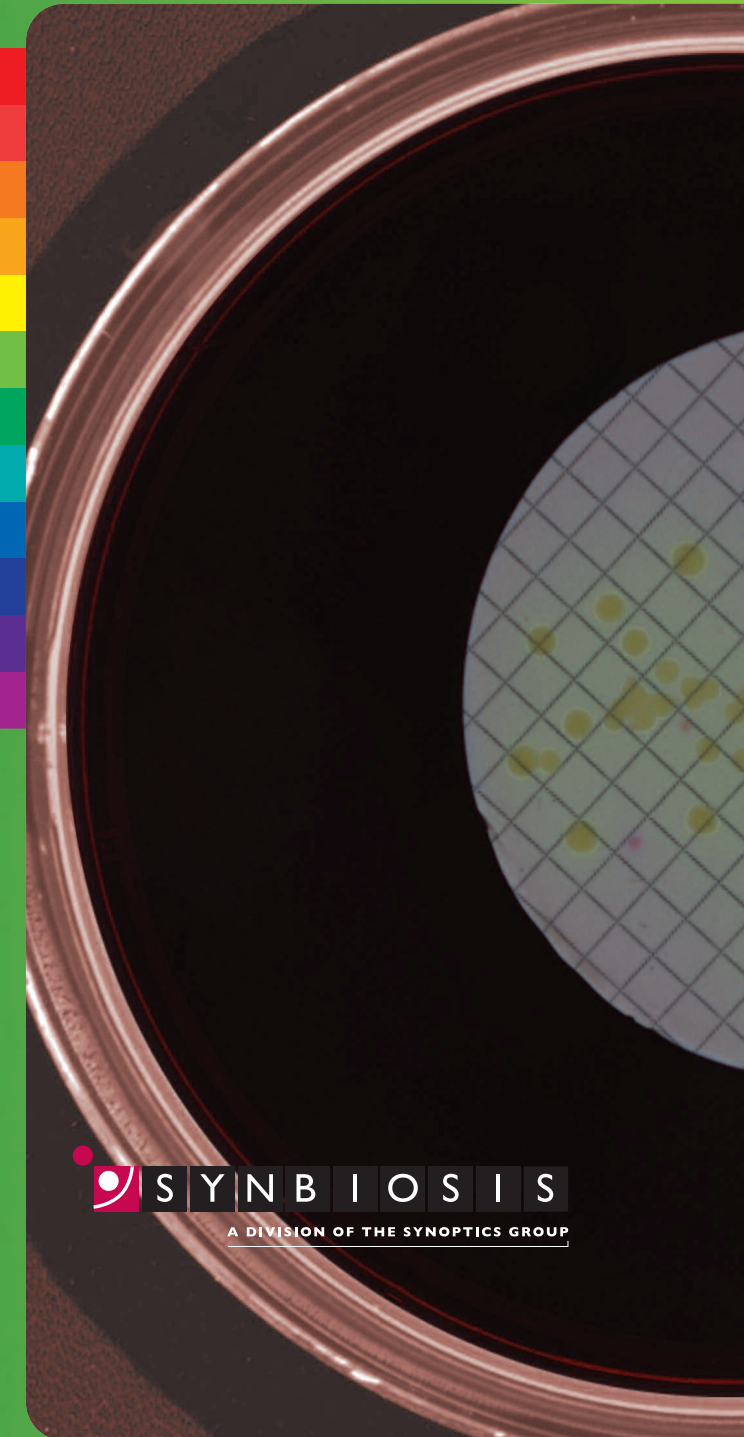

A DIVISION OF THE SYNOPTICS GROUP

MICROBIAL LIMITS TEST-MEMBRANE FILTRATION

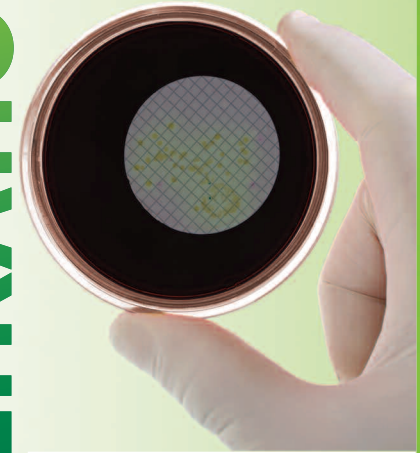

COUNTS ON GRIDDED MEMBRANE



DILUTION FACTOR



COUNTS FOR EACH COLONY COLOUR



REPORTING



ProtoCOL 2 with its easy to use software, built in MS SQL server database and touch screen interface makes the counting of Membrane filters fast, secure and accurate

- Microbial Limits Test is designed to perform the quantitative estimations of specific viable microorganisms present in samples
- There are four methods for this test, including membrane filtration
- Membrane filtration applies to samples which contain antimicrobial substances
- Samples are filtered through a membrane filter which retains any microorganisms present. The filters are placed on a suitable growth medium and incubated. Any viable aerobic organisms will grow and be counted