

ProtoCOL 2

Colony counting and zone sizing system

Quick Start Guide - connecting the touch screen monitor

ProtoCOL 2 does not have any antivirus software installed. Add own antivirus software if required

1. Unpack the instrument and the monitor and place on a secure level surface.
2. Attach the black bracket to the instrument via the silver mounting plate as shown below. The four screws are secured using a Philips screw driver (not supplied).



3. Remove the protective wrapping from the cables.
4. With the monitor facing towards the front of the instrument connect the cables as shown below.



5. Slide the bracket that is attached to the monitor over the top of the bracket that is now attached to the front of Protocol 2 taking care to feed the cables under the monitor so as not to impede the movement of the monitor.

-
6. Secure the monitor onto the bracket using the 4 screws.



CONTACT SYNBIOSIS:

EUROPE
BEACON HOUSE
NUFFIELD ROAD
CAMBRIDGE
CB4 1TF

Tel: +44 (0)1223 727125

Fax: +44 (0)1223 727101

Email: eurosales@synbiosis.com

Email: intlsales@synbiosis.com

www.synbiosis.com

IN USA:

5108 PEGASUS COURT, SUITE M
FREDERICK
MD 21704

Tel: 800 686 4451 (toll free) / 301 662 2863

Fax: 301 631 3977

Email: ussales@synbiosis.com