


ProtoCOL 2- Installing Honeywell Voyager Scanner 9520-38-3

1 Unpacking the scanner

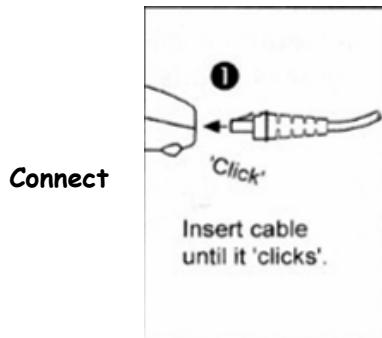

➔ Unpack the scanner and connect the 10-pin RJ45 male connector of the USB cable into the jack on the scanner. There will be an audible *click* when the connection is made


2 Installing barcode software



Run

➔ Run the install Barcode exe

3 Connecting the scanner

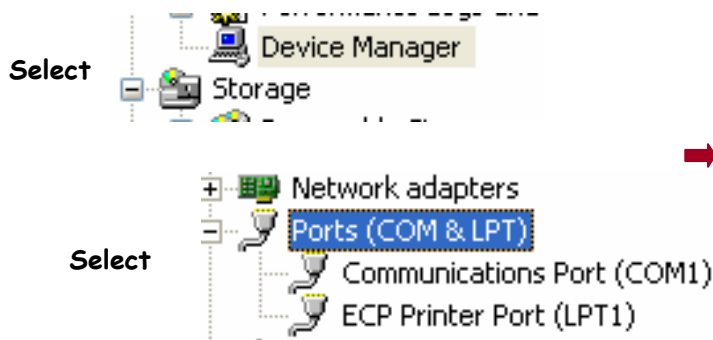

➔ Connect the scanner to a free USB port


(Optional) Click on the Control Panel, Administrative Tools, Computer Management and open the Device Manager and Ports (COM & LPT), the scanner will **not** yet be visible

4 Scan the barcode



➔ Scan the barcode to enable bi-directional USB emulation

5 Device Manager

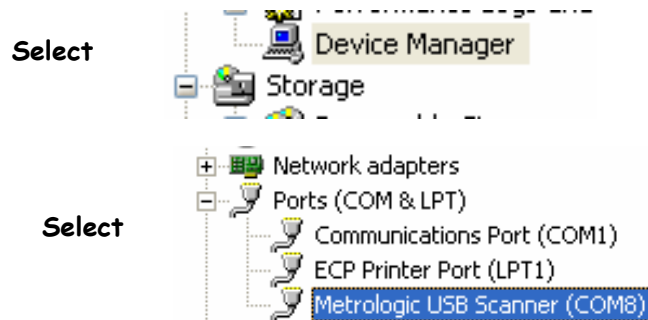

➔ (Optional) Open Windows Device Manager- *Ports (COM & LPT)*, the scanner will now appear as a Metrologic USB Scanner. *Note - the COM port allocation will vary*

6 Select barcode



After completing colour classification, click the Measure tab and select Use Barcodes. Accept the batch

The scanned plate barcode will appear in the plate ID box

ProtoCOL 2 is now ready to make the first measurement

CONTACT SYNBIOSIS:
EUROPE:
Tel: +44 (0)1223 727125

USA:
Tel: 800 686 4451/301 662 2863

Email: support@synbiosis.com

Email: support@synbiosis.com

www.synbiosis.com
P2HVS1210