

ProtoCOL 2- Editing a plate count - Quick Guide

1 Measuring a plate

Click




➔ After measuring the plate click on the line of results that require editing. The Edit Result tab will now appear

2 Editing a plate

Click



➔ Click on the Edit Result tab and the edit functions will appear

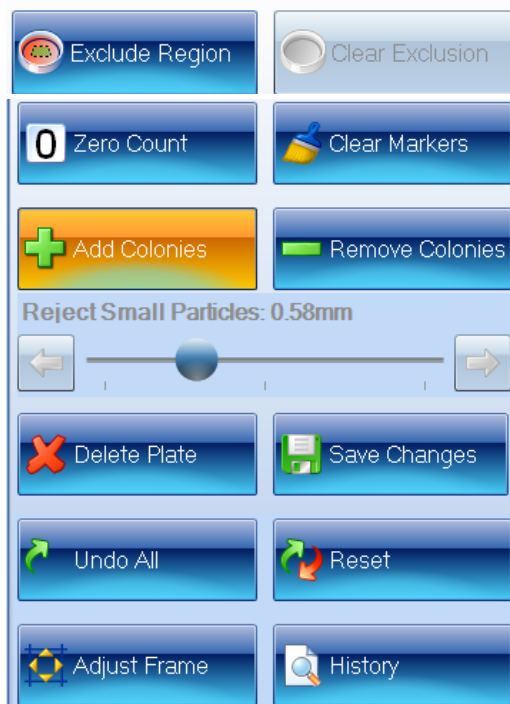

➔ Edit the plate count as required using the edit functions


Click



➔ Click Save Changes


3 Enter a reason for the edit

Select

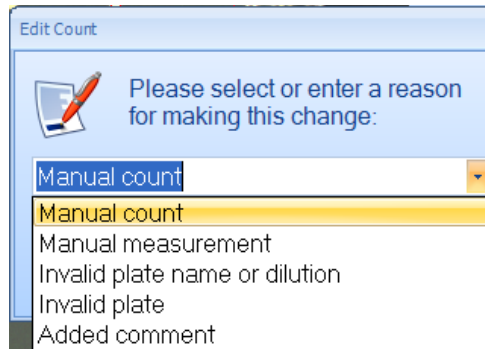

➔ Add a reason or select from the dropdown list the reason for making the change

Click

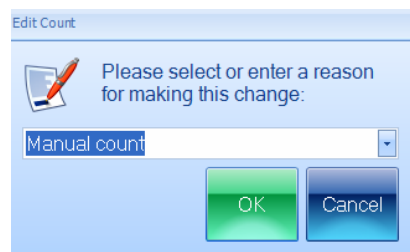

➔ Click OK

4 Results

Plate Name	Diluti...	Colony Name	Count / Frame	Coun...	Mean Co...	User	Flags	Created	Comments /...
123	1:1	A	102	122	102	P2ADMIN	M	04/04/2011 11:18:28	

➔ The edited results will now appear in the results table with a flag to show an adjustment has been made

CONTACT SYNBIOSIS:
EUROPE:
Tel: +44 (0)1223 727125
Email: support@synbiosis.com

USA:
Tel: 800 686 4451/301 662 2863
Email: support@synbiosis.com

www.synbiosis.com
P2EPC0411