

ProcScan - Setting Up Quick Guide

1 Install scanner software

- ➔ Install the scanner software onto ProtoCOL 2 or PC according to scanner instructions
- ➔ Connect USB cable to ProtoCOL 2 or PC

2 Start ProtoCOL 2 software



- ➔ Log onto ProtoCOL 2 software

3 Twain source

Click




- ➔ Click on the ProtoCOL 2 icon in the top left of the screen

Select



- ➔ From the drop down menu select Settings then "Configure External TWAIN Source"

Select

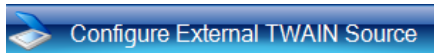

- ➔ Select the scanner being used


4 Select scanner

Click



- ➔ Select the Image tab

Click

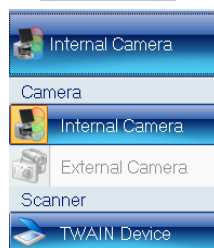

- ➔ Select the scanner by selecting TWAIN Device from the drop down menu on the left side of the screen labelled Internal Camera if scanner is being used with a ProtoCOL 2

5 Capture image

Click



- ➔ Click Capture Image
- ➔ The scanner software will now open and the settings can be set for the type of plate to be analysed

CONTACT SYNBIOSIS:
EUROPE:
Tel: +44 (0)1223 727125
Email: support@synbiosis.com

USA:
Tel: 800 686 4451/301 662 2863
Email: support@synbiosis.com

www.synbiosis.com
PSSUQ61210