

ProtoCOL 2 spiral plates - setting up a dilution Quick Guide

1 Create a spiral batch

Create



→ Create and accept a spiral batch on ProtoCOL 2

2 Enter dilution

Click



→ Click on the dilution button

Select




→ Select Edit Dilution List

3 Enter dilution series

Click



→ Click Add

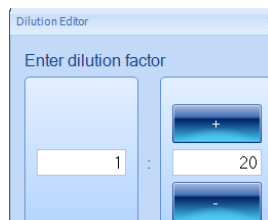

→ Enter dilution by either clicking on +/- and dilutions increase by a factor of 10. Alternatively you can click in text box and enter dilution

Click



→ After each dilution is added click Add

→ If a dilution needs removing, highlight dilution requiring deletion and click Remove

Click



→ When the dilution series is complete click Close

4 Select dilution



→ The dilutions can now be selected from the drop-down menu or by scrolling through using the arrow keys. When the required dilution is selected it will be displayed in the dilution box and the measurement can be taken

ProtoCOL 2 is now ready to make the first measurement

CONTACT SYNBIOSIS:
EUROPE:
Tel: +44 (0)1223 727125
Email: support@synbiosis.com

USA:
Tel: 800 686 4451/301 662 2863
Email: support@synbiosis.com

www.synbiosis.com
P2SPSUD0111