

ProtoCOL 3 - Exporting/Importing Batches Quick Guide

1 Create a batch

➔ Create a batch as per relevant quick guide

2 Export batch

Click




➔ Click on ProtoCOL 3 icon

Select



➔ Select Export Data as XML

Click

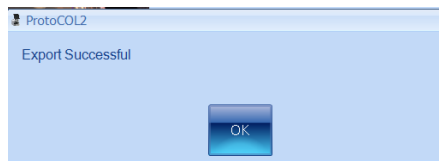

➔ Set an export path and name the export batch

➔ Click OK

3 Create a batch based on exported batch

Select




➔ Click on New Batch

Select



➔ Click on Create from Exported Batch

Click

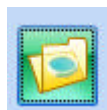

➔ Click on the file icon to locate the exported batch

Please select a batch file to import:

test export.xml



Click




➔ Click Go

4 Test measure



➔ Click Test Measure


5 Enter a plate ID

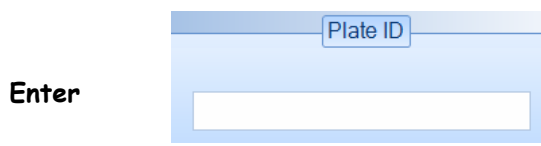


➔ Enter a plate ID found on Measure tab. Auto-increment default is "on", to change this click on the Configuration tab



6 Name the batch

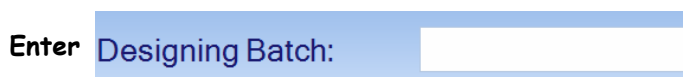

7 Accept batch

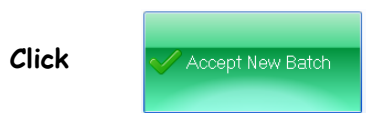

➔ Accept batch

ProtoCOL 3 is now ready to make the first measurement

CONTACT SYNBIOSIS:
EUROPE:
Tel: +44 (0)1223 727125
Email: support@synbiosis.com

USA:
Tel: 800 686 4451/301 662 2863
Email: support@synbiosis.com

www.synbiosis.com
P3EIB1211