


ChromaZona

MIC Strip Batch Set Up - Quick Guide

1

Start ChromaZona software

Enter

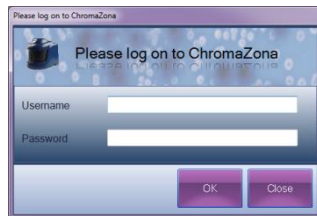

➔ Log on to ChromaZona

2

Position plate

➔ Insert the plate holder, ensuring the plate holder which is recommended by the software is used

➔ Place plate onto plate holder

3

Create batch

Click



➔ Click New Batch

4

Plate configuration

Click

➔ Defaults are Circular plate 90mm
Change if necessary

5

Choose batch type

Select



➔ Select MIC Strip

6

Name the Batch

Enter



7

Select the Measure Tab

Select



8

Enter a Plate ID

Enter




Enter a plate ID



Auto-increment default is “on”, to change this click on the Configuration tab

9

Save the batch

Click



Click the Accept New Batch button to save the selected settings

10

Select the MIC Strip Supplier


Select



11

Detect Strips

Click

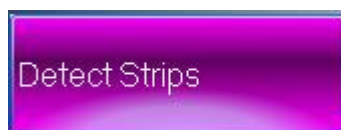

Select the Detect Strips button to automatically detect the MIC strips present on the agar plate

12

Measure Plate

Click



CONTACT SYNBIOSIS:

EUROPE:
Tel: +44 (0)1223 727125
Email: support@synbiosis.com

USA:
Tel: 800 686 4451/301 662 2863
Email: support@synbiosis.com