

# ChromaZona

## Software Update - Quick Guide

### 1 Start PC


➔ Log onto PC as an Administrator and shut all running software down

### 2 Insert software source

➔ Insert the software source, unzipping if the installation set is zipped. Run the file Setup.exe and follow the on-screen instructions

3 The PC may require a Windows update (particularly if the PC is Windows XP, Windows Vista or Windows 7). If you are presented with a message stating “[This application needs Microsoft .Net Framework 4.0 full version which is not yet installed. This can be found in the Support directory.](#)”, then follow step 3 if you have a full installation set or 4 if you just have an update set

#### Support folder

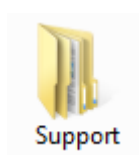

➔ Open the Support folder



➔ Open the Framework folder



➔ Run SetupFramework.exe. Follow the on-screen instructions, performing a system restart if update required

➔ Run Setup.exe once again and follow the on-screen instructions (see ChromaZona Full Installation Guide for details)

4

## Download



dotNetFx40\_Full\_x86\_x64.exe

Open a web browser, download and run the .NET Framework 4.0 full update from:

<http://www.microsoft.com/en-gb/download/details.aspx?id=17718>

5



Setup.exe

Run Setup.exe once again and follow the on-screen instructions (see ChromaZona Full Installation Guide for details)

**ChromaZona is now ready to be used**

### CONTACT SYNBIOSIS:

EUROPE:

Tel: +44 (0)1223 727125

Email: [support@synbiosis.com](mailto:support@synbiosis.com)

USA:

Tel: 800 686 4451/301 662 2863

Email: [support@synbiosis.com](mailto:support@synbiosis.com)

[www.synbiosis.com](http://www.synbiosis.com)

CZSU0716