


ProtoCOL 3

Entering Plate IDs by CVS - Quick Guide

1 Enter a plate list

Create

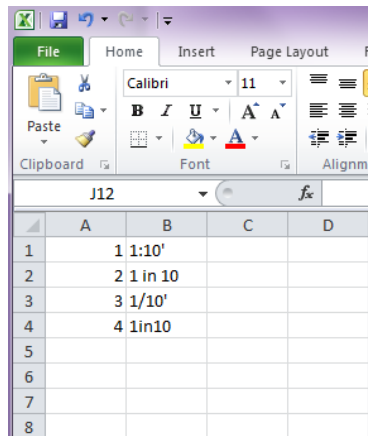

	A	B	C	D
1	1	1:10'		
2	2	1 in 10		
3	3	1/10'		
4	4	1in10		
5				
6				
7				
8				

Create a plate list in Excel and save it as a CSV file. The plate list needs to be in columns and can include plate IDs in one column and dilutions in quotation marks in a second column. The dilutions can be in any of the formats shown

2 Start ProtoCOL 3 software

Enter



Please log on to ProtoCOL3

Please log on to ProtoCOL3

Username

Password

OK Close

Log on to ProtoCOL using the username P3Admin and the password is the ProtoCOL 3 serial number

3 Create batch

Click



Create a new batch, after completing the classification click on the configuration tab


Select



Select Configuration tab

4

Import plate list

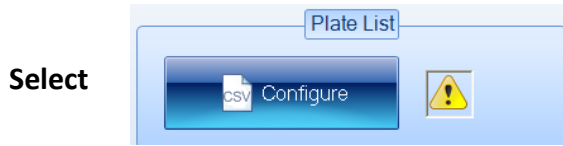

Select Configure


Select Import CSV File from the source for plate IDs

Click Open button to select CSV file to import

Verify plate list, if satisfactory click OK

5

Accept batch

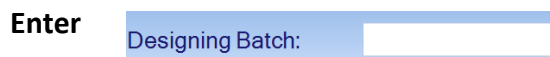

Name the Batch



Accept batch

ProtoCOL 3 is now ready to measure the first plate in the plate list

CONTACT SYNBIOISIS:

EUROPE:
Tel: +44 (0)1223 727125
Email: support@synbiosis.com
www.synbiosis.com

USA:
Tel: 800 686 4451/301 662 2863
Email: support@synbiosis.com
P3IDSCVS1216