

ProtoCOL 3

Exporting/Importing Batches - Quick Guide

1

Create a batch

➔ Create a batch as per relevant quick guide

Export batch

Click



➔ Click on ProtoCOL 3 icon


2

Select



➔ Select Export Data as XML

Click

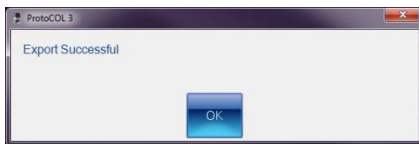

➔ Set an export path and name the export batch

➔ Click OK

Create a batch based on exported batch

Select



➔ Click on New Batch

Select




➔ Click on Create from Exported Batch

3

Click



➔ Click on the file icon to locate the exported batch

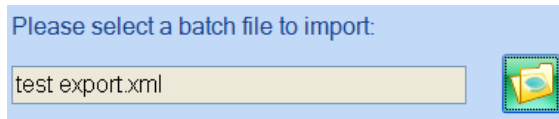

Click



➔ Click Go

Test measure

4

Click



Click Test Measure

Enter a plate ID

5

Enter




Enter a plate ID found on Measure tab. Auto-increment default is "on", to change this click on the Configuration tab

Click



Name the batch

6

Enter




Name the batch

Accept batch

7

Click



Accept batch

ProtoCOL 3 is now ready to make the first measurement

CONTACT SYNBIOSIS:

EUROPE:
Tel: +44 (0)1223 727125
Email: support@synbiosis.com
www.synbiosis.com

USA:
Tel: 800 686 4451/301 662 2863
Email: support@synbiosis.com
P3EIB1216