


# ProtoCOL 3

## Printing Plate Layouts - Quick Guide

### 1 Layouts

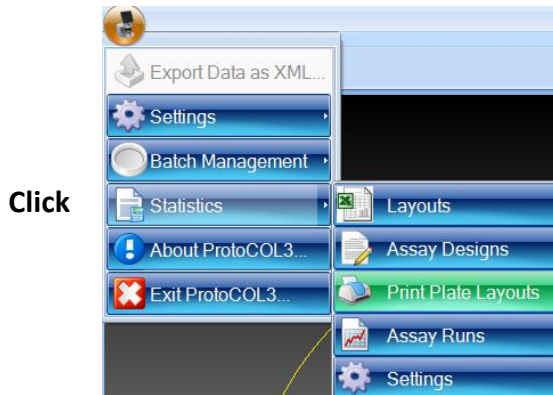


Click

Click on the ProtoCOL 3 icon in top left of screen

Select Statistics

Select Print Plate Layouts

### 2 Select the Layout to Print

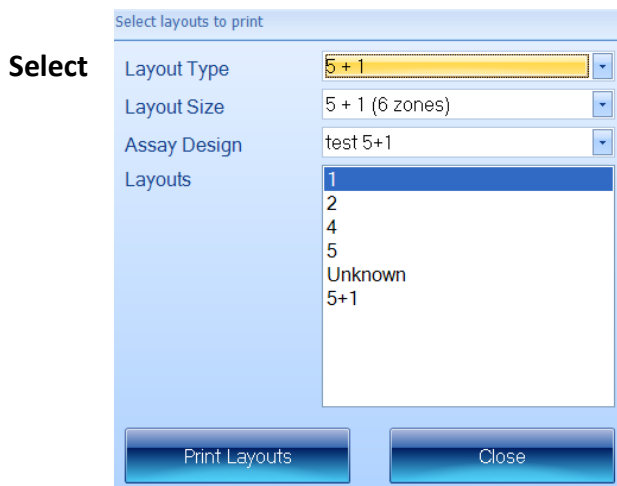

Select

Select Layout Type and Size from dropdown

Select Assay Design from dropdown

Highlight the Layout for printing

Click Print Layouts

Click Close

**CONTACT SYNBIOISIS:**

EUROPE:  
 Tel: +44 (0)1223 727125  
 Email: [support@synbiosis.com](mailto:support@synbiosis.com)  
[www.synbiosis.com](http://www.synbiosis.com)

USA:  
 Tel: 800 686 4451/301 662 2863  
 Email: [support@synbiosis.com](mailto:support@synbiosis.com)  
**P3PPL1216**