

ProtoCOL 3

Re-measuring a plate from a plate list


This Quick Guide explains how to delete a measured plate from a pre-defined plate list and allow that plate name to be used again in the measurement step

1

Plate Name	Diluti...	Colony Name	Count / Frame
4	1 : 1	A	244
3	1 : 1	A	60
2	1 : 1	A	261
1	1 : 1	A	262


➔ In the 'Results' tab click on the plate you want to re-measure and go to the 'Edit/Review' tab

2

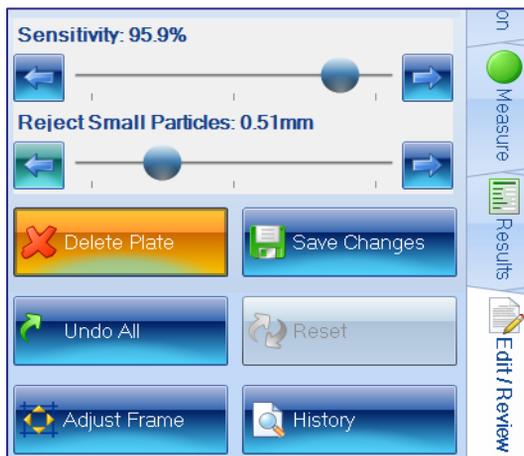

➔ Select 'Delete Plate' then 'Save Changes'

3

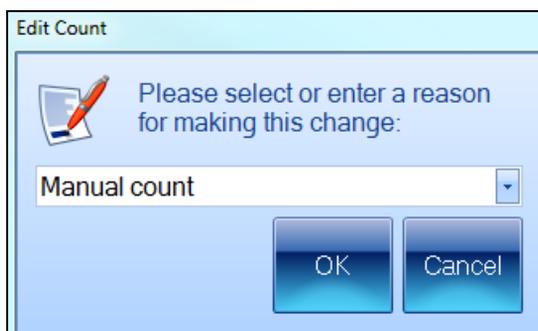


➔ Choose an audit reason from the dropdown menu or enter your own

4




➔ A dialog box will open asking if you would like to be able to re-use that plate name for measuring - select 'Yes' if you want to re-measure this plate

5

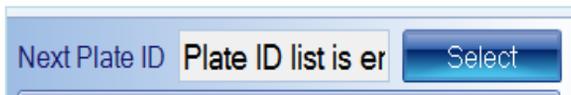

➔ Press 'Select' in the 'Next Plate ID' box to view the plate names that are available to measure.

6

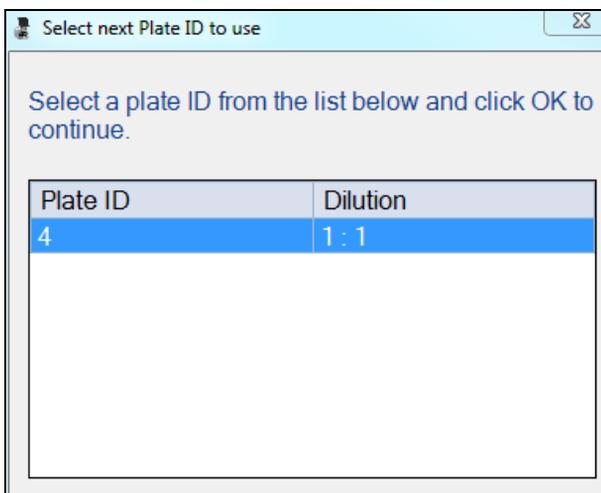

➔ Choose the plate you want to re-measure from the plate list and press OK

You can now re-measure this plate

CONTACT SYNBIOISIS:

EUROPE:
 Tel: +44 (0)1223 727125
 Email: support@synbiosis.com
www.synbiosis.com

USA:
 Tel: 800 686 4451/301 662 2863
 Email: support@synbiosis.com
P3RMP0818