

ProtoCOL 3

Connecting Touch Screen PC - Quick Guide

1 Unpack ProtoCOL 3

➔ Unpack the instrument and the PC and place on a secure level surface

Attach bracket

2



➔ Attach the black bracket to the instrument via the silver mounting plate as shown below. The four screws are secured using a pozidrive screwdriver (not supplied)

Attach monitor

3



- ➔ Remove the protective wrapping from the cables
- ➔ With the PC facing towards the front of the instrument connect the cables as shown below
- ➔ Slide the bracket that is attached to the monitor over the top of the bracket that is now attached to the front of ProtoCOL taking care to feed the cables under the monitor so as not to impede the movement of the monitor
- ➔ Secure the monitor onto the bracket using the 4 screws

THE LOGIN DETAILS FOR THE PROTOCOL 3 PLUS ARE
Username: P3ADMIN
Password: The unit's serial number in capitals

CONTACT SYNBIOSIS:

EUROPE:
Tel: +44 (0)1223 727125
Email: support@synbiosis.com
www.synbiosis.com

USA:
Tel: 800 686 4451/301 662 2863
Email: support@synbiosis.com
P3CTSPC1216