

Protos 3 - Chromogenic ID - Batch set-up Quick Guide

1 Start Protos 3 software

Click



➔ Run Protos 3

2 Position plate

➔ Insert the plate holder, ensuring the plate holder which is recommended by the software is used

➔ Place plate onto plate holder

3 Create batch

Click



➔ Click New Batch

4 Plate configuration

Click



➔ Defaults are Circular plate 90mm
Change if necessary

5 Choose batch type

Select



➔ Select Chromogenic Media

6 Enter a Plate ID

Enter



➔ Enter a plate ID found on Measure tab. Autoincrement default is "on", to change this click on the Configuration tab

Click



7

Save the batch

Enter

Designing Batch:



Name the Batch

Click



Accept batch

Protos 3 is now ready to make the first measurement

CONTACT SYNBIOSIS:
EUROPE:
Email: support@synbiosis.com
www.synbiosis.com
Protos3CBQG