

aCOLyte software installation - Quick Guide

1 Log onto PC

- ➔ Log onto PC as Administrator
- ➔ Plug the aCOLyte into a mains socket


2 Run software installation



- ➔ Run the software installation
- ➔ Click Next to start installation process
- ➔ The Device Driver Installation Wizard dialog box will be displayed, click Next
- ➔ When the device driver installation is complete press Finish


3

Destination

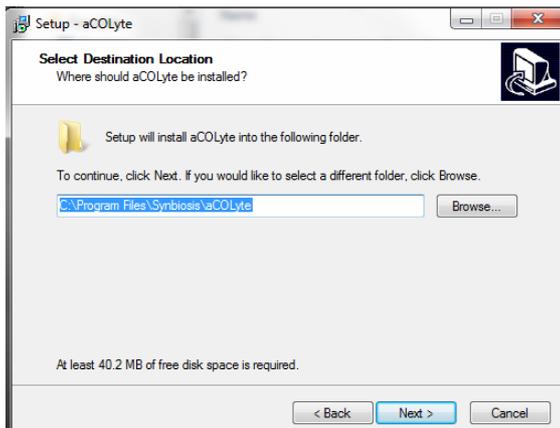

➔ Select a destination where aCOLyte should be installed

➔ Select Next

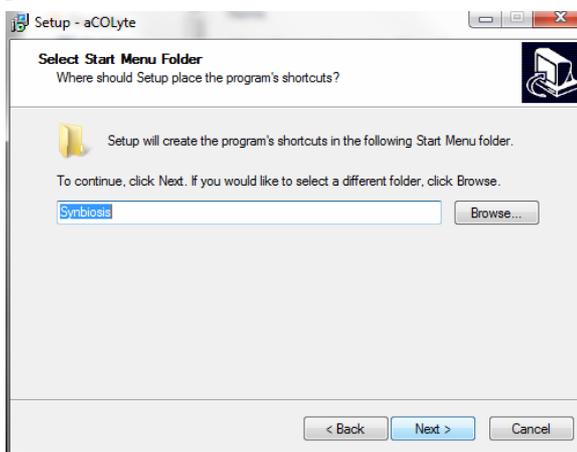


➔ Select a place for program shortcuts

➔ Select Next

4

Installation

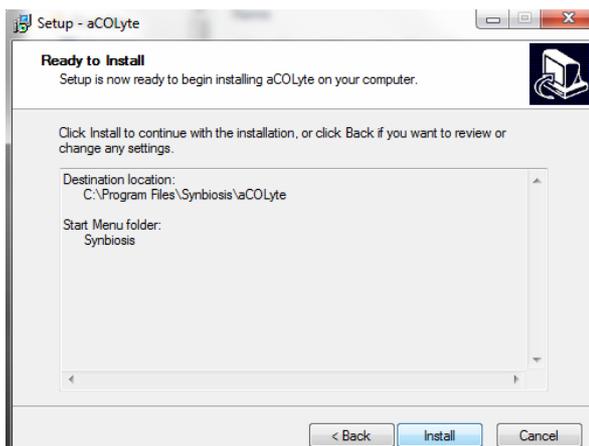


➔ Select Install

➔ Select Finish


5

Security Code

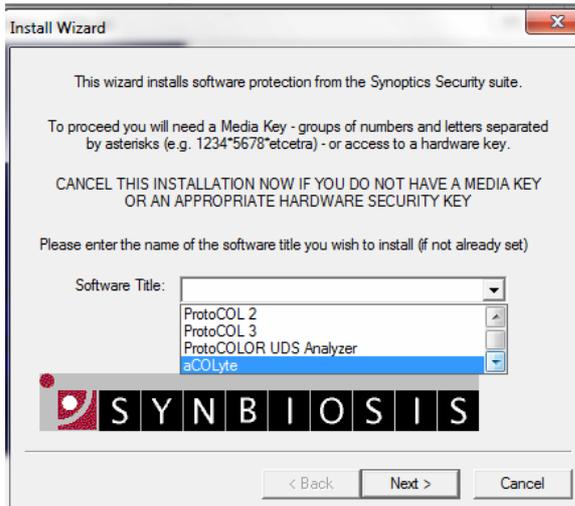


➔ Select aCOLyte

➔ Click Next

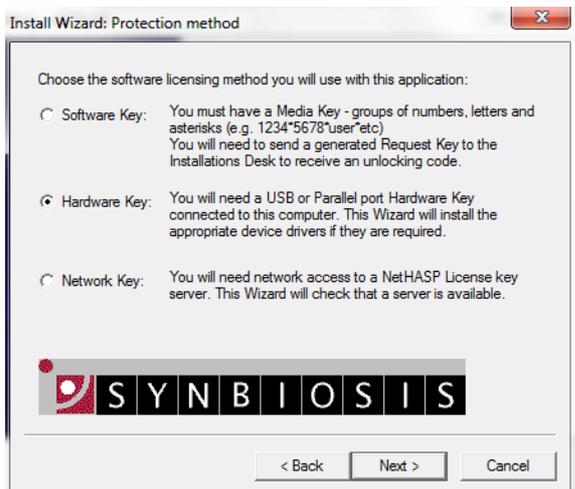

➔ Select the key appropriate to your system


➔ Click Next

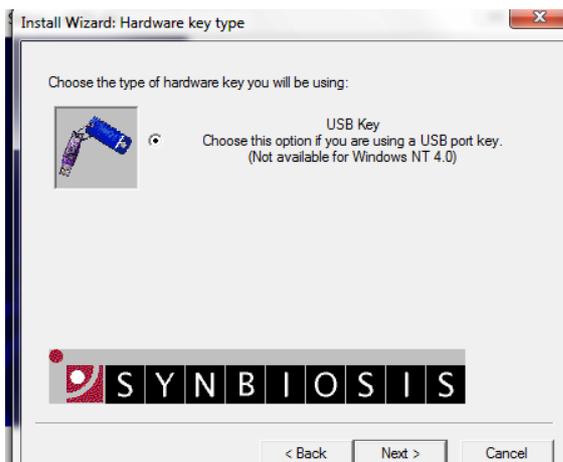

➔ If a hardware key is being used, select the type of hardware that will be used

➔ Click Next

➔ Prepare the PC to install the device driver for the security key. Make sure there are no security keys attached. Click Next to install drivers - this may take a few seconds

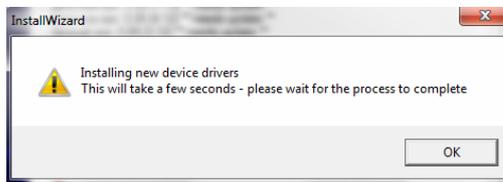

➔ Click OK



➔ Connect the hardware key and follow the instructions from the dialogue box

➔ Click Next



➔ When you are told that the test was successful click Finish to exit or Back to install the security key on other USB PORT

➔ Plug in the USB lead and double click the aCOLyte; the software should now open.

CONTACT SYNBIOSIS:
EUROPE:
Tel: +44 (0)1223 727125
Email: support@synbiosis.com

USA:
Tel: 800 686 4451/301 662 2863
Email: support@synbiosis.com

www.synbiosis.com
ACOLSI0612



A DIVISION OF THE SYNOPTICS GROUP